

NOTE NO. 17 - January 15 1987

THERAPEUTIC RANGE DETERMINATION OF 3,4-DIAMINOPYRIDINE

Analytes: 3,4-Diaminopyridine (100 ng/ml)


Sample Matrix: Human serum

Sample Size: 10 microliters

Column: 5 micron GFF ISRP, 5 cm x 4.6 mm ID

Regis Product Number: 731450

Mobile Phase: 98% 0.01 M KH_2PO_4 (pH 6.8), 2% Methanol
(HPLC grade reagents)

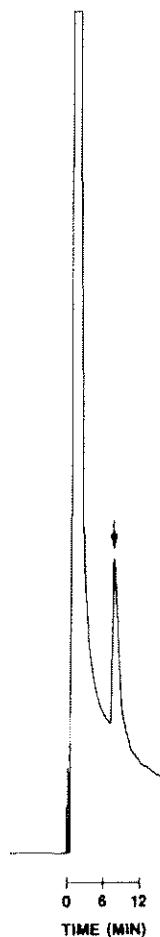

Detection: 295 nm, 0.002 AUFS

Flow Rate: 1.0 ml/min

Note: Sample filtered through 0.2 um membrane prior to direct injection

T.J. Szczerba, Regis Chemical Company,
8210 Austin Avenue, Morton Grove, IL 60053