

NOTE NO. 10 - July 1 1986

THERAPEUTIC RANGE DETERMINATION OF PROPRANOLOL

Analytes: Propranolol (100 ng/ml)

Sample Matrix: Human serum

Sample Size: 10 microliters

Column: 5 micron GFF ISRP, 5 cm x 4.6 mm ID

Regis Product Number: 731450

Mobile Phase: 90% 0.1 M KH_2PO_4 (pH 6.8), 10 % Acetonitrile
(HPLC grade reagents)

Detection: 290 nm, 0.001 AUFS

Flow Rate: 2.0 ml/min

Note: Sample filtered through 0.2 μm
membrane prior to direct injection



**T.J. Szczerba and J.A. Perry, Regis Chemical Company,
8210 Austin Avenue, Morton Grove, IL 60053**